



Tutela V4 - Release Notes

<u>Version</u>	<u>Description</u>	<u>Release Date</u>
01	Issue	15/06/2016

Contents

- 1.0 Overview 2
- 2.0 New Features 2
- 4.0 Removed/Retired Features 3
- 5.0 Bug Fixes 3
- 6.0 System Requirements 3

1.0 Overview

1.1 This 4th generation website has been developed in line with the principles of Tutela Version 3.2 website which is specifically designed to meet the requirements of the highly regulated Medical and Pharmaceutical markets for the validation, management and activity audit tracking of electronic data records.

1.2 This version as with all previous versions provides remote access to the automatic recorded data and alarm exceptions of the sensing probes installed in the customer's facility. Customers can gain access to their data via the secure Tutela V4 web base application hosted by Tutela medical. The data records are published in real time via these secure web pages. Remote site hosting, maintenance and validation of electronic data records are regulated through Tutela Medical **UK ISO 9001 Business Management System**.

2.0 New Features

- 2.1 Downloading of raw data in Excel format.
- 2.2 Automatic identification of country during login.
- 2.3 Password reset functionality.
- 2.4 Pin code reset functionality.
- 2.5 Automatically move probes to the top of the sensor list with outstanding incidents.
- 2.6 User management system.
- 2.7 Contact management system.
- 2.8 Access to view and manage probe isolations.
- 2.9 Manage alarm contacts and lists.
- 2.10 Digital Contact types integrated into system. (configurable feature)
- 2.11 Display a fault instead of a value, if a probe has not recorded data for a set period.
- 2.12 Auto scaling of graphs.
- 2.13 Incident report summary of individual probes.
- 2.14 Graphs to link to incident audit.
- 2.15 Site audits can be viewed by website function and search with calendar functionality.
- 2.16 Export Isolation History.
- 2.17 Validate password length to 8 characters strong.

3.0 Improvements/Enhancements

- 3.1 New Style look and feel.
- 3.2 More flexible date range.
- 3.3 New PDF library functionality.
- 3.4 Toggle between alarm and alert settings pages.

4.0 Removed/Retired Features

- 4.1 Validate icon removed from sensor list page.

5.0 Bug Fixes

- 5.1 Dynamic Daylight saving to remove issue of viewing data in the current daylight saving zone only.

6.0 System Requirements

- 6.1 Operating Systems and browsers compatibility is as follows:

Yes
No

	Windows 10	Windows 7 64-Bit	Windows 7	Windows XP
Edge		N/A	N/A	N/A
IE 11			N/A	N/A
IE 10	N/A		N/A	N/A
IE 9 x 64	N/A			N/A
IE 9	N/A			N/A
IE 8	N/A			
Chrome 51				
Chrome 50				
Chrome 49				
Chrome 48		N/A		
Chrome 47		N/A		
Chrome 46		N/A		
Chrome 45		N/A		
Chrome 44		N/A		
Chrome 43		N/A		
Chrome 42	N/A	N/A		

Chrome 41	N/A	N/A		
FireFox 47				
FireFox 46				
FireFox 45				
Firefox 44	N/A	N/A		
Firefox 43	N/A	N/A		
Firefox 42	N/A	N/A		
Firefox 41	N/A	N/A	N/A	N/A
Firefox 40	N/A	N/A	N/A	N/A
Firefox 39	N/A	N/A	N/A	N/A
Firefox 38	N/A	N/A	N/A	N/A
Opera 37			N/A	N/A
Opera 36				
Opera 35				
Safari 9	N/A	N/A	N/A	N/A
Safari 8	N/A	N/A	N/A	N/A
Safari 7	N/A	N/A	N/A	N/A
Safari 6.2	N/A	N/A	N/A	N/A

	MAC 10.11	MAC 10.10	MAC 10.9	MAC 10.8
Edge	N/A	N/A	N/A	N/A
IE 11	N/A	N/A	N/A	N/A
IE 10	N/A	N/A	N/A	N/A
IE 9 x 64	N/A	N/A	N/A	N/A
IE 9	N/A	N/A	N/A	N/A
IE 8	N/A	N/A	N/A	N/A
Chrome 51				
Chrome 50				
Chrome 49				
Chrome 48				
Chrome 47				
Chrome 46				
Chrome 45				
Chrome 44				
Chrome 43				
Chrome 42				
Chrome 41				
FireFox 47				
FireFox 46				
FireFox 45				
Firefox 44				
Firefox 43				
Firefox 42				



Firefox 41	N/A	N/A	N/A	N/A
Firefox 40	N/A	N/A	N/A	N/A
Firefox 39	N/A	N/A	N/A	N/A
Firefox 38	N/A	N/A	N/A	N/A
Opera 37				N/A
Opera 36				
Opera 35				
Safari 9			N/A	N/A
Safari 8	N/A		N/A	N/A
Safari 7	N/A	N/A		N/A
Safari 6.2	N/A	N/A	N/A	