

Connected Automated Monitoring + v4.2.2
(Formerly known as Tutela)
Release Notes

CONFIDENTIALITY

The information contained within this manual is proprietary. Ownership is solely vested in Checkit UK Limited.

This information may not be copied or reproduced or passed on to any third party without the prior written permission of Checkit UK Limited.

The Checkit Tutela system Hardware and Software whether whole or in part are subject to the laws of copyright.

The adoption and/or use of techniques functionality and know how embodied in this document, the hardware and/or software source and/or runtime code modules may not be copied or utilized in any way by third parties without the prior written consent of Checkit UK Limited.

<u>Version</u>	<u>Description</u>	<u>Release Date</u>
01	Issue	15/06/2016
1.1	Issue	29/09/2016
1.2	Issue	03/02/2017
1.3	Issue	04/07/2017
1.4	Issue	03/11/2017
1.5	Issue	01/03/2018
1.6	Issue	10/10/2018
2.0	Issue	01/06/2020
2.1	Issue	21/09/2020
2.2	Issue	09/11/2020

Contents

1.0 Overview	5
2.0 Additional Features.....	5
3.0 Improvements/Enhancements	5
4.0 Removed/Retired Features.....	7
5.0 Bug Fixes	7
6.0 System Requirements	7

1.0 Overview

1.1 This document outlines the additional features, Improvements, and bug fixes to the 4th generation website.

1.2 This version as with all previous versions provides remote access to the automatic recorded data and alarm exceptions of the sensing probes installed in the customer's facility. Customers can gain access to their data via the secure Connected Automated Monitoring web base application. The data records are published in real time via these secure web pages. Remote site hosting, maintenance and validation of electronic data records are regulated through Tutela Medical UK ISO 9001 Business Management System.

2.0 Additional Features


3.0 Improvements/Enhancements

1. Setpoint values can no longer be set outside the specified limits. An error message will be shown if tried to do so.

The screenshot displays the 'Connected Automated Monitoring +' web interface. The header includes the Checkit logo, the user 'imran.khurram@checkit.net', and navigation links for 'SITE LIST', 'EXIT', 'REPORT', 'HELP', and 'LOGOUT'. The main content area shows 'ALARM SETTINGS' for 'NEXTCSL - Next Control Systems - IT Monitoring'. A table lists channels with their descriptions, current values, and alarm setpoints. A modal dialog titled 'Change setpoint' is open, showing a text input field with '70.0 °C' and a message: 'Please enter a new value between 0.0 and 35.0'. The dialog has 'CONFIRM' and 'EXIT' buttons. A red box in the top right corner highlights an information icon and the message 'Value is outside the limits!'.

CHANNEL	DESCRIPTION	VALUES	LOW ALARM SETPOINT	DELAY	HIGH ALARM SETPOINT	DELAY
Control Chart	Testing	42.9 °C	0.0 °C	0.0 min	0.0 °C	0.0 min
Server Ambient	Server Room Ambient	21.4 °C	0.0 °C	0.0 min	12.0 °C	0.0 min

2. User will be locked if he/she tried to login with wrong password three times.

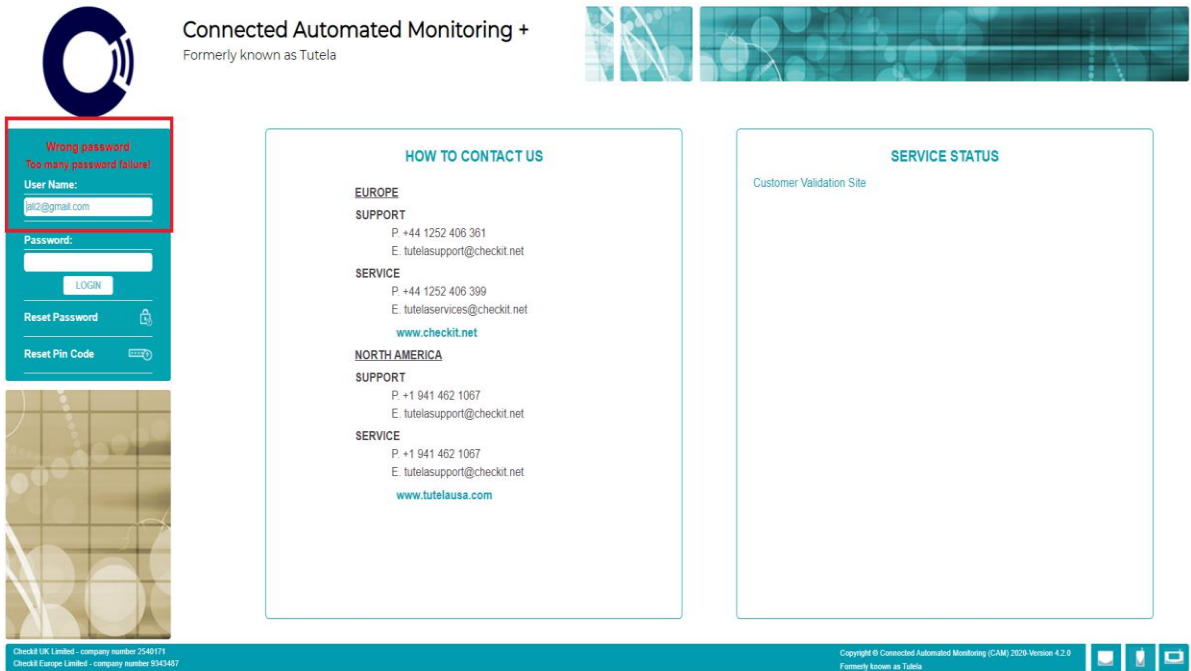

Wrong password
Too many password failures!

User Name:
ali2@gmail.com

Password:
[input field]

LOGIN

Reset Password

Reset Pin Code

HOW TO CONTACT US

EUROPE

SUPPORT
P: +44 1252 406 361
E: tutelasupport@checkit.net

SERVICE
P: +44 1252 406 399
E: tutelaservices@checkit.net
www.checkit.net

NORTH AMERICA

SUPPORT
P: +1 941 462 1067
E: tutelasupport@checkit.net

SERVICE
P: +1 941 462 1067
E: tutelasupport@checkit.net
www.tutelausa.com


SERVICE STATUS

Customer Validation Site

Checkit UK Limited - company number 2540171
Checkit Europe Limited - company number 9343407

Copyright © Connected Automated Monitoring (CAM) 2020-Version 4.2.0
Formerly known as Tutela

This locked user can now be enabled by Admins from “System User Status” page. Upon enabling the user, the wrong password failure count will be reset, and user can login again.

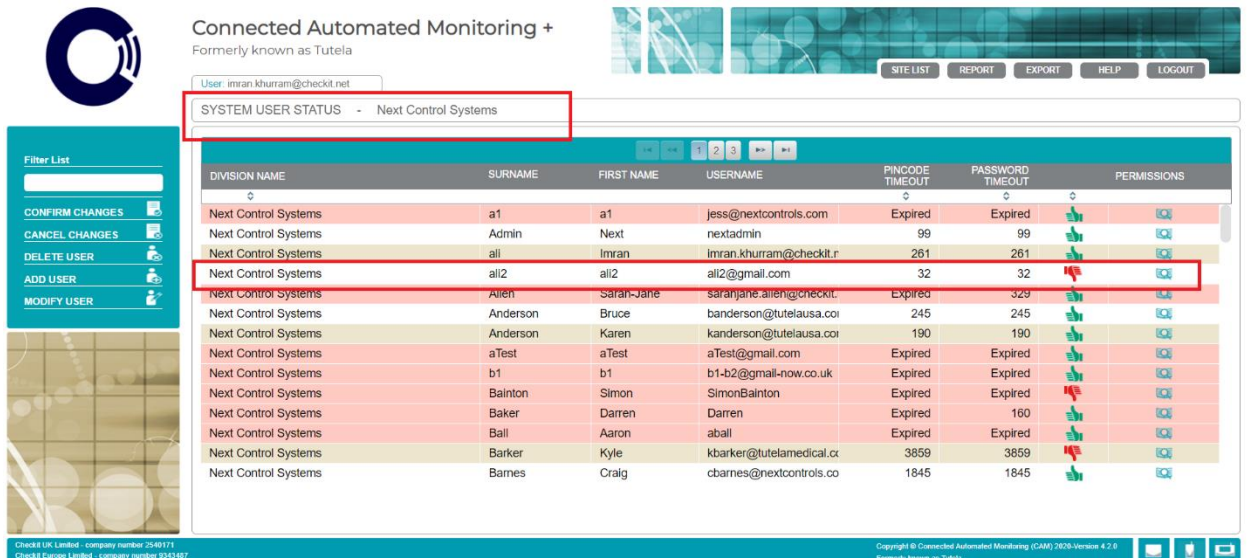

Connected Automated Monitoring +
Formerly known as Tutela

User: imran.khurram@checkit.net

SYSTEM USER STATUS - Next Control Systems

DIVISION NAME	SURNAME	FIRST NAME	USERNAME	PINCODE TIMEOUT	PASSWORD TIMEOUT	PERMISSIONS
Next Control Systems	a1	a1	jess@nextcontrols.com	Expired	Expired	[lock icon]
Next Control Systems	Admin	Next	nextadmin	99	99	[lock icon]
Next Control Systems	ali	Imran	imran.khurram@checkit.r	261	261	[lock icon]
Next Control Systems	ali2	ali2	ali2@gmail.com	32	32	[lock icon]
Next Control Systems	Alien	Sarah-Jane	sarahjane.aliens@checkit.	Expired	329	[lock icon]
Next Control Systems	Anderson	Bruce	banderson@tutelausa.co	245	245	[lock icon]
Next Control Systems	Anderson	Karen	kanderson@tutelausa.co	190	190	[lock icon]
Next Control Systems	aTest	aTest	aTest@gmail.com	Expired	Expired	[lock icon]
Next Control Systems	b1	b1	b1-b2@gmail-now.co.uk	Expired	Expired	[lock icon]
Next Control Systems	Bainton	Simon	SimonBainton	Expired	Expired	[lock icon]
Next Control Systems	Baker	Darren	Darren	Expired	160	[lock icon]
Next Control Systems	Ball	Aaron	aball	Expired	Expired	[lock icon]
Next Control Systems	Barker	Kyle	kbarker@tutelamedical.co	3859	3859	[lock icon]
Next Control Systems	Barnes	Craig	cbarnes@nextcontrols.co	1845	1845	[lock icon]

Checkit UK Limited - company number 2540171
Checkit Europe Limited - company number 9343407

Copyright © Connected Automated Monitoring (CAM) 2020-Version 4.2.0
Formerly known as Tutela

4.0 Removed/Retired Features

5.0 Bug Fixes

6.0 System Requirements

Yes
No

	Windows 10	Windows 8.1	Windows 8	Windows 7 64-Bit	Windows 7	Windows XP
Edge 81		N/A	N/A	N/A	N/A	N/A
Edge 80		N/A	N/A	N/A	N/A	N/A
Edge 79		N/A	N/A	N/A	N/A	N/A
Edge 18		N/A	N/A	N/A	N/A	N/A
Edge 17		N/A	N/A	N/A	N/A	N/A
Edge 16		N/A	N/A	N/A	N/A	N/A
Edge 15		N/A	N/A	N/A	N/A	N/A
IE 11			N/A		N/A	N/A
IE 10	N/A	N/A			N/A	N/A
IE 9 x64	N/A	N/A	N/A		N/A	N/A
IE 9	N/A	N/A	N/A			N/A
IE 8	N/A	N/A	N/A	N/A		
Chrome 81					N/A	N/A
Chrome 80					N/A	N/A
Chrome 79					N/A	N/A
Chrome 78					N/A	N/A
Chrome 77					N/A	N/A
Chrome 76					N/A	N/A
Chrome 75					N/A	N/A
Chrome 74					N/A	N/A
Chrome 73					N/A	N/A
Chrome 72						N/A
Chrome 71						N/A
Chrome 70						N/A
Chrome 69						N/A
Chrome 68						N/A
Chrome 67						N/A
Chrome 66						N/A
Chrome 65						N/A



Chrome 64						N/A
Chrome 63						N/A
Chrome 62						N/A
Chrome 61						N/A
Chrome 60						N/A
Chrome 59						N/A
Chrome 58						N/A
Chrome 57						N/A
Chrome 56						N/A
Chrome 55						N/A
Chrome 54						N/A
Chrome 53						N/A
Chrome 52						N/A
Chrome 51						N/A
Chrome 50						N/A
Chrome 49						
Chrome 48						
Chrome 47						
Chrome 46						
Chrome 45						
Chrome 44						
Chrome 43						
Chrome 42	N/A					
Chrome 41	N/A					
FireFox 76		N/A	N/A		N/A	N/A
FireFox 75			N/A		N/A	N/A
FireFox 74		N/A	N/A		N/A	N/A
FireFox 73		N/A	N/A		N/A	N/A
FireFox 72					N/A	N/A
FireFox 71					N/A	N/A
FireFox 70					N/A	N/A
FireFox 69					N/A	N/A
FireFox 68					N/A	N/A
FireFox 67					N/A	N/A
FireFox 66					N/A	N/A
FireFox 65						N/A
FireFox 64						N/A
FireFox 63						N/A
FireFox 62						N/A
FireFox 61						N/A
FireFox 60						N/A
FireFox 59						N/A
FireFox 58						N/A



Firefox 57						N/A
Firefox 56						N/A
Firefox 55						N/A
Firefox 54						N/A
Firefox 53						N/A
Firefox 52						N/A
Firefox 51						N/A
Firefox 50						N/A
Firefox 49						N/A
Firefox 48						N/A
Firefox 47						N/A
Firefox 46						
Firefox 45						
Firefox 44						
Firefox 43						
Firefox 42						
Firefox 41						
Firefox 40						
Firefox 39						
Firefox 38						
Opera 68					N/A	N/A
Opera 67					N/A	N/A
Opera 66					N/A	N/A
Opera 65					N/A	N/A
Opera 64					N/A	N/A
Opera 63					N/A	N/A
Opera 60					N/A	N/A
Opera 58					N/A	N/A
Opera 57					N/A	N/A
Opera 56					N/A	N/A
Opera 55					N/A	N/A
Opera 54						N/A
Opera 53					N/A	N/A
Opera 52					N/A	N/A
Opera 51					N/A	N/A
Opera 50					N/A	N/A
Opera 49					N/A	N/A
Opera 48						N/A
Opera 47						N/A
Opera 46						N/A
Opera 45						N/A
Opera 44						N/A
Opera 43						N/A
Opera 42						N/A



Opera 41						N/A
Opera 40						N/A
Opera 39						N/A
Opera 38						N/A
Opera 37					N/A	N/A
Opera 36						
Opera 35						
Safari 13	N/A	N/A	N/A	N/A	N/A	N/A
Safari 12	N/A	N/A	N/A	N/A	N/A	N/A
Safari 11	N/A	N/A	N/A	N/A	N/A	N/A
Safari 10.1	N/A	N/A	N/A	N/A	N/A	N/A
Safari 9	N/A	N/A	N/A	N/A	N/A	N/A
Safari 8	N/A	N/A	N/A	N/A	N/A	N/A
Safari 7	N/A	N/A	N/A	N/A	N/A	N/A
Safari 6.2	N/A	N/A	N/A	N/A	N/A	N/A

	Mac 10.15	Mac 10.14	Mac 10.13	Mac 10.12	Mac 10.11	Mac 10.10	Mac 10.9	Mac 10.8
Edge 80		N/A	N/A	N/A	N/A	N/A	N/A	N/A
Edge 79		N/A	N/A	N/A	N/A	N/A	N/A	N/A
Chrome 81					N/A	N/A	N/A	N/A
Chrome 80 x64						N/A	N/A	N/A
Chrome 79						N/A	N/A	N/A
Chrome 78 x64						N/A	N/A	N/A
Chrome 77 x64	N/A						N/A	N/A
Chrome 76 x64	N/A						N/A	N/A
Chrome 75 x64	N/A						N/A	N/A
Chrome 74 x64	N/A						N/A	N/A
Chrome 73 x64	N/A						N/A	N/A
Chrome 72 x64	N/A						N/A	N/A
Chrome 71 x64	N/A						N/A	N/A
Chrome 70 x64	N/A						N/A	N/A
Chrome 69 x64	N/A						N/A	N/A
Chrome 68 x64	N/A	N/A					N/A	N/A
Chrome 67 x64	N/A	N/A					N/A	N/A
Chrome 66 x64	N/A	N/A					N/A	N/A
Chrome 65 x64	N/A	N/A					N/A	N/A
Chrome 64 x64	N/A	N/A					N/A	N/A
Chrome 63 x64	N/A	N/A					N/A	N/A
Chrome 62 x64	N/A	N/A					N/A	N/A
Chrome 61 x64	N/A	N/A					N/A	N/A



Chrome 60 x64	N/A	N/A					N/A	N/A
Chrome 59 x64	N/A	N/A					N/A	N/A
Chrome 58 x64	N/A	N/A					N/A	N/A
Chrome 57 x64	N/A	N/A					N/A	N/A
Chrome 56 x64	N/A	N/A					N/A	N/A
Chrome 55 x64	N/A	N/A					N/A	N/A
Chrome 54 x64	N/A	N/A	N/A				N/A	N/A
Chrome 53 x64	N/A	N/A	N/A				N/A	N/A
Chrome 52 x64	N/A	N/A	N/A				N/A	N/A
Chrome 51 x64	N/A	N/A	N/A	N/A			N/A	N/A
Chrome 50 x64	N/A	N/A	N/A	N/A			N/A	N/A
Chrome 49 x64	N/A	N/A	N/A	N/A				
Chrome 48 x64	N/A	N/A	N/A	N/A				
Chrome 47 x64	N/A	N/A	N/A	N/A				
Chrome 46 x64	N/A	N/A	N/A	N/A				
Chrome 45 x64	N/A	N/A	N/A	N/A				
Chrome 44 x64	N/A	N/A	N/A	N/A				
Chrome 43 x64	N/A	N/A	N/A	N/A				
Chrome 42 x64	N/A	N/A	N/A	N/A				
Chrome 41 x64	N/A	N/A	N/A	N/A				
FireFox 76		N/A	N/A	N/A	N/A	N/A	N/A	N/A
FireFox 75					N/A	N/A	N/A	N/A
FireFox 74					N/A	N/A	N/A	N/A
FireFox 73						N/A	N/A	N/A
FireFox 72					N/A	N/A	N/A	N/A
FireFox 71 x64					N/A	N/A	N/A	N/A
FireFox 70						N/A	N/A	N/A
FireFox 69	N/A						N/A	N/A
FireFox 68	N/A						N/A	N/A
FireFox 67	N/A						N/A	N/A
FireFox 66	N/A						N/A	N/A
FireFox 65	N/A						N/A	N/A
FireFox 64	N/A						N/A	N/A
FireFox 63	N/A						N/A	N/A
FireFox 62	N/A						N/A	N/A
FireFox 61	N/A	N/A					N/A	N/A
FireFox 60	N/A	N/A					N/A	N/A
FireFox 59	N/A	N/A						N/A
FireFox 58	N/A	N/A					N/A	N/A
FireFox 57	N/A	N/A						N/A
FireFox 56	N/A	N/A						N/A
FireFox 55	N/A	N/A	N/A					N/A
FireFox 54	N/A	N/A	N/A					N/A
FireFox 53	N/A	N/A	N/A					N/A



FireFox 52	N/A	N/A	N/A					N/A
FireFox 51	N/A	N/A	N/A					N/A
FireFox 50	N/A	N/A	N/A					N/A
FireFox 49	N/A	N/A	N/A					N/A
FireFox 48	N/A	N/A	N/A					
FireFox 47	N/A	N/A	N/A					
FireFox 46	N/A	N/A	N/A	N/A				
FireFox 45	N/A	N/A	N/A	N/A				
FireFox 44	N/A	N/A	N/A	N/A				
FireFox 43	N/A	N/A	N/A	N/A				
FireFox 42	N/A	N/A	N/A	N/A				
FireFox 41	N/A	N/A	N/A	N/A				
FireFox 40	N/A	N/A	N/A	N/A				
FireFox 38	N/A	N/A	N/A	N/A				
Opera 66		N/A				N/A	N/A	N/A
Opera 65		N/A				N/A	N/A	N/A
Opera 64						N/A	N/A	N/A
Opera 63	N/A						N/A	N/A
Opera 62	N/A						N/A	N/A
Opera 60	N/A						N/A	N/A
Opera 58	N/A						N/A	N/A
Opera 57	N/A						N/A	N/A
Opera 56	N/A						N/A	N/A
Opera 55	N/A	N/A					N/A	N/A
Opera 54	N/A	N/A					N/A	N/A
Opera 53	N/A	N/A					N/A	N/A
Opera 52	N/A	N/A					N/A	N/A
Opera 51	N/A	N/A					N/A	N/A
Opera 50	N/A	N/A					N/A	N/A
Opera 49	N/A	N/A						N/A
Opera 48	N/A	N/A						N/A
Opera 47	N/A	N/A	N/A					N/A
Opera 46	N/A	N/A	N/A				N/A	N/A
Opera 45	N/A	N/A	N/A					N/A
Opera 44	N/A	N/A	N/A					N/A
Opera 43	N/A	N/A	N/A					N/A
Opera 42	N/A	N/A	N/A					N/A
Opera 41	N/A	N/A	N/A					N/A
Opera 40	N/A	N/A	N/A					N/A
Opera 39	N/A	N/A	N/A					N/A
Opera 38	N/A	N/A	N/A					N/A
Opera 37	N/A	N/A	N/A	N/A				N/A
Opera 36	N/A	N/A	N/A	N/A				
Opera 35	N/A	N/A	N/A	N/A				



Safari 13		N/A	N/A	N/A	N/A	N/A	N/A	N/A
Safari 12	N/A		N/A	N/A	N/A	N/A	N/A	N/A
Safari 11	N/A	N/A		N/A	N/A	N/A	N/A	N/A
Safari 10	N/A	N/A	N/A		N/A	N/A	N/A	N/A
Safari 9	N/A	N/A	N/A	N/A		N/A	N/A	N/A
Safari 8	N/A	N/A	N/A	N/A	N/A		N/A	N/A
Safari 7	N/A	N/A	N/A	N/A	N/A	N/A		N/A
Safari 6.2	N/A	N/A	N/A	N/A	N/A	N/A	N/A	